Wasser-Destillierautomaten

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Automatic water distillation apparatus for the production of distilled water, corresponding to the European Pharmaceutical Code, with a conductivity below $2,3 \mu\text{S}/\text{cm}$ at $+20^\circ\text{C}$, depending on the condition of the raw water.

By actuating the main switch all operating courses required for the distillation process are controlled fully automatically, i.e. water supply, regulation of cooling water, switching on of heating, switching off of heating in case of lack of water, switching off of the heating and the cooling water supply when the storage vessel is filled, and switching on when distillate has been taken. The respective operating position is indicated by signal lamps.

By utilization of the cooling water for distillate production heated up at 80°C , heating energy and water are saved.

Thermostatic low water cut-off protects the heating element in case of water shortage.

Condenser (cooling coil) interchangeable.

The apparatus are made of special "stainless steel" 1.4301, Heating Element are made of "stainless steel" No 1.4876.

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Elektroanschluß 230V~, 50/60 Hz, Anschlußwert 0,22kVA.

Decalcifying pump RO 1060

for the removal of water sediments in automatic water distillation apparatus. The cleaning of the still, of the condenser and of the piping system is done in one operating cycle by pumping decalcifying solution in circulation through the system.

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RO 221065

Decalcifying agent

for removing lime and magnesium deposits. Packing unit 5 kg.

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Für Destillierautomaten
RO 1032-1034

Separate supply unit RO 1044

for the separate feed of the still with completely demineralized water. As no water heated up in the condenser is utilized, the output of the automatic distillation apparatus is reduced by approx. 15%.

Water pressure: min. 0.2 bar, max. 10 bar, electrical connection 230V~, 50/60 Hz.

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RO 553 111

Phosphate cartridge RO 1051

which treats the cooling water, prevents the recrystallization of the hardness causing salts. Applicable for all automatic distillation apparatus which are operated with separate supply 1044. For installation into the water supply pipe.

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Dechlorit filter RO 1053

removes the chlorine added by the waterworks. Suitable for use with all distillation apparatus, for installation in the water inlet pipe.

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Ec declaration of conformity according to the CE Tension Directive and EMC Directive.

Herewith we declare that ROWA Water Bath models:
R0 03003, R0 03042, R0 03043, R0 03044

Pressureless Steam Sterilizer model:
R0 04120

with the following technical data:

Nominal voltage 230 a.c., 50 / 60 Hz

Protection class I

Protection type covered

The apparatus comply with the requirements of the German Equipment Safety and were tested according to:

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ROWA Labortechnische Anlagen GmbH